


VDAS E-LAB STUDENT QUICKSTART GUIDE

VERSATILE DATA ACQUISITION SYSTEM E-LAB




Collect data from a remotely performed experiment


1. If it is not already installed, download the TecEquipment VDAS@e-lab software from:
<https://www.tecequipment.com/downloads>
2. Extract the VdasSetup_xxx.exe from the downloaded zip file and double click it to run.
3. Agree to the licensing conditions and click 'Install'

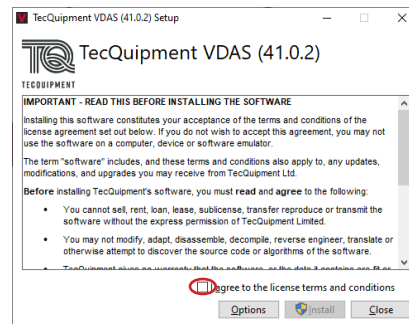


4. Wait for the Installation to complete, then select Launch.

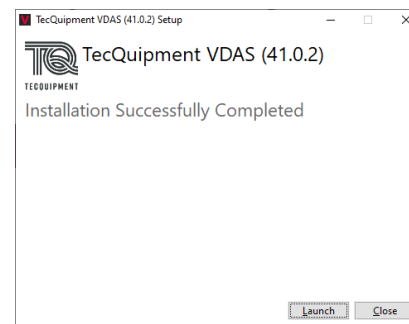

5. If the software has already been installed:
Launch it by clicking on the new desktop icon or start menu item:

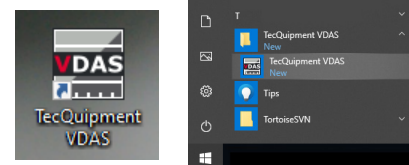


6. Select Options, Select VDAS Application:
7. Select the name of the apparatus being used:

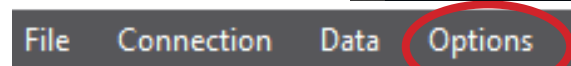


8. Select Connection, Communication Settings




9. Select the Remote Participant tab:
10. Then click on: Connect to a shared experiment:

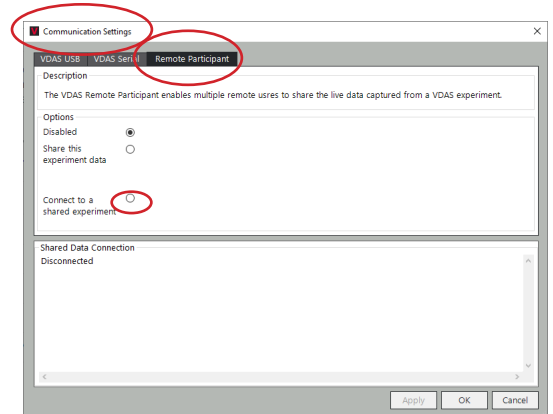

11. Enter your Shared ID (provided by your tutor).
12. Enter your name (this is your registration as being present for the experiment).
13. Once correctly connected the Status box will contain a tick.
14. When connected click OK to return to the data window.

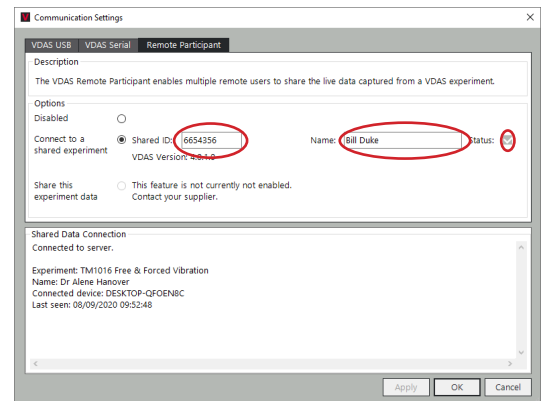

15. Experiment data is now being shared. You can independently start viewing, recording and analysing data via the VDAS tool bar controls.

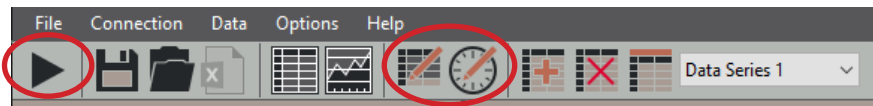

Access VDAS online help at any time by pressing F1.